



**NANO**●**PARTZ**<sup>™</sup>

The gold nanoparticle for nanotechnology

# **Conjugating Carboxyl- Polymer Spherical Gold Nanoparticles to Antibodies Kit**

Part Number 21-PC-KIT1 Rev 1.1  
Product Profile

©Nanopartz, All Rights Reserved  
3/1/2009

## Purpose

Conjugation, purification, and validation of conjugating Carboxyl-polymer SGNPs to Goat anti-Human.

## Introduction

This is a kit to instruct the conjugation of carboxyl-polymer spherical gold nanoparticles (SGNPs) to any primary amine using EDC chemistry. This kit includes IgG Goat anti-Human as the conjugation target.

EDC is a carboxyl and amine-reactive zero-length crosslinker. EDC reacts with a carboxyl group first and forms an amine reactive *O*-acylisourea intermediate that quickly reacts with an amino group to form an amide bond and release of an isourea by-product (see the Additional Information Section). The intermediate is unstable in aqueous solutions, and therefore, two step conjugation procedures require *N*-hydroxysuccinimide for stabilization.<sup>1,2</sup> Failure to react with an amine will result in hydrolysis of the intermediate, regeneration of the carboxyl and release of an *N*-substituted urea. A side reaction is the formation of an *N*-acylurea, which is usually restricted to carboxyls located in hydrophobic regions of proteins.<sup>1,3</sup>

## Kit Applications

- Conjugate carboxyl to amine groups in peptides, proteins, and DNA labeling through 5' phosphate groups

## Kit Includes

1. 20-PC-KIT1-1 Carboxyl-polymer SGNPs to IgG Conjugation Kit Instruction Manual
2. 20-PC-KIT1-2 20-PC-xx-50, Conjugated SGNPs in MES, Carboxyl-polymer, xx nm, 50 OD mLs, where xx is the size of the nanoparticle delivered
3. 20-PC-KIT1-3 *N*-ethyl-*N'*-dimethylaminopropyl-carbodiimide (EDC), 10 mg in 1.5 ml microcentrifuge tube
4. 20-PC-KIT1-4 Nanopartz PBS based Buffer, pH=7.4, 10 ml
5. 20-PC-KIT1-5 IgG Goat Anti-Human, 10 µl (1 mg/ml solution)
6. 20-PC-KIT1-6 18 MEG DI H<sub>2</sub>O, 10 ml
7. 20-PC-KIT1-7 Low binding 1.5 ml microcentrifuge tubes

## Storage and Shelf Life

Refrigerate at 4°C. Shelf life is 14 days, limited by the IgG. Use only the EDC amount necessary – once dissolved in water its lifetime is less than one hour. Allow kit components to warm to room temperature prior to using.

# NANOPARTZ™

The gold nanoparticle for nanotechnology

## **Other required materials/instruments**

1. Transfer pipettes for volumes ranging from 1  $\mu$ l to 2 ml
2. Microcentrifuge with minimum rcf = 10,000 for 20 nm sizes and larger, for smaller sizes, please contact us for the rcf necessary
3. Vortex for microcentrifuge tubes
4. Sonicator

## **Other optional materials/instruments**

1. Malvern Nano ZS DLS
2. UV VIS

## Supplementary Information

EDC reacts with a carboxyl group first and forms an amine-reactive *O*-acylisourea intermediate that quickly reacts with an amino group to form an amide bond and release of an isourea by-product (Figure 1).

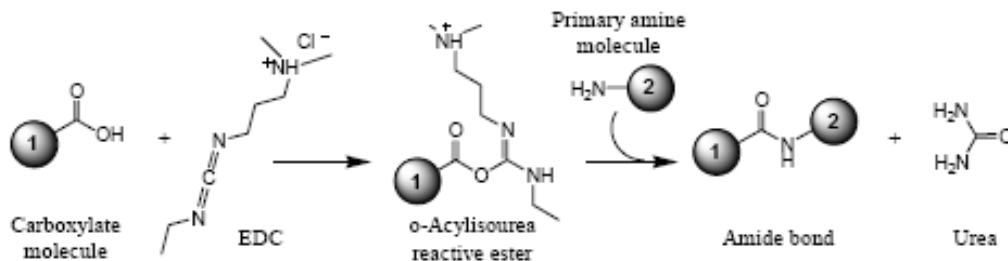


Figure 1 One-step EDC reaction with carboxyl and amine-containing molecules

## References

1. Grabarek, Z. and Gergely, J. (1990). Zero-length crosslinking procedure with the use of active esters. *Anal Biochem* 185:131-5.
2. Staros, J.V., *et al.* (1986). Enhancement by *N*-hydroxysulfosuccinimide of water-soluble carbodiimide-mediated coupling reactions. *Anal Biochem* 156:220-2.
3. Timkovich, R. (1977). Detection of the stable addition of carbodiimide to proteins. *Anal Biochem* 79:135-43.

This product ("Product") is warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation") and to be free from defects in material and workmanship. Unless otherwise expressly authorized in writing, Products are supplied for research use only. No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to one year from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than the original purchaser of the Product ("Buyer").

**No other warranties, express or implied, are granted, including without limitation, implied warranties of merchantability, fitness for any particular purpose, or non infringement. Buyer's exclusive remedy for non-conforming Products during the warranty period is limited to replacement of or refund for the non-conforming Product(s).**

There is no obligation to replace Products as the result of (i) accident, disaster or event of force majeure, (ii) misuse, fault or negligence of or by Buyer, (iii) use of the Products in a manner for which they were not designed, or (iv) improper storage and handling of the Products.

© Nanopartz, All Rights Reserved 2008

Nanopartz, A Division of Concurrent Analytical Inc.  
www.nanopartz.com